

## Release Notes

Proteograph™ Analysis Suite v2.0  
August 3, 2022

- **PAS Proteogenomics Workflow to identify and explore peptides with protein sequence altering genetic variants**
  - **Build a custom peptide database**
    - New feature to build a customized peptide sequence database using user-uploaded custom or sample-specific variant call files (VCF).
      - VCF files are used to annotate genetic variants that may result in amino acid variants (i.e., variant peptides) not captured in the canonical reference database.
    - Variant peptide sequences are combined with the reference database to generate customized peptide sequence databases.
  - **Search for variant peptides**
    - Using the customized peptide sequence database, now users can search their DDA data for variant peptides using MSFragger.
  - **Variant Peptide Browser and Proteogenomics Data Explorer**
    - *Variant Peptide Browser* – Identified variant peptides are summarized in interactive plots and table.
      - Summary table shows all identified variant peptides and can be sorted/filtered by sample, nanoparticle, protein, variant, and allele frequency. Each variant (row) links to a view within the Proteogenomic Data Explorer.
      - The number of variant peptides found across samples, conditions, and nanoparticles are shown. The distribution of variant peptide intensity is stratified by reference and alternative alleles and homozygous/heterozygous alleles across samples, conditions, and nanoparticles.
    - *Proteogenomics Data Explorer* – Interactive tool to explore peptide and variant peptide data maps in genomic space for entire proteins and at nucleic acid/amino acid resolution
      - Gene structure, protein structure, domain information, and functional region information are displayed.
      - Amino acid variants within variant peptides are highlighted, including both reference and alternative alleles.
- **Improvements to analysis protocols**
  - Integration of MSFragger database search engine.
    - Pre-installed MSFragger-based Proteogenomics protocol (DDA, VCF required).
  - Enabled DIA-NN MBR feature<sup>1</sup>.
- **Dashboard to monitor PAS account usage**
  - Track the total number of plates added, projects added, completed analyses, and data storage.
  - Added links to most recently added plates, projects, completed analyses, and data storage.
- **Expansion and improvement of analysis visualizations**
  - Added additional background datasets and overlay for multiple samples in waterfall plots.
  - Added ability to interact with sample correlation and similarity plots and visualize plots showing underlying data for each comparison.
  - Added option to view and compare all metrics from plate map grid across all wells and across all plates analyzed.

- Added option to select order of comparison for differential abundance group analysis.
- Added new interactivity with differential abundance group analysis volcano plot display.
- **Changes to file and project management**
  - Added feature to multi-select projects when launching analyses.
  - Added feature to collapse Analyses folders by project.
  - Added project filtering capabilities.
- **Improvements to analysis stability and plot visualizations for large-scale study sizes**
  - Stability and browser performance upgrades.
  - Analysis visualization and plot rendering speed improvements.
  - Streamlined analysis summary menu improves navigation between visualizations.
- **Added PAS software end user license agreement (EULA)**
- **Changes to help content**
  - Updated user guide.
  - Updated tooltips.
- **Other general improvements**
  - Sample Description File is now stored in Data Files.
  - Minor bug fixes.

1. Demichev, V., *et al.* DIA-NN: Neural Networks and interference correction enable deep proteome coverage in high throughput. *Nature Methods* **17**, 41–44 (2019).

© 2022 Seer, Inc. Seer, Proteograph, and the Seer logo are trademarks of Seer, Inc. MSFragger © 2016 The Regents of the University of Michigan. All other marks are the property of their owners. For Research Use Only. Not for use in diagnostic procedures. Use of the Proteograph Analysis Suite is subject to the terms and conditions contained in Seer's end user license agreement. Use of PAS may incur separate cloud compute and data storage fees.